

Fluorometer Quick Installation Guide

- 1. Unpack:** Open the packaging box, which contains the instrument, packing list, and accessories as shown in Figure 1.

NOTE: *The instrument should be installed on a clean, dry, and level surface for proper operation.*

- 2. Product Inspection:** Please inspect the components of the product according to the items listed in the "Packing List," as shown in Figure 2, including the instrument, power cord, power adapter, USB drive (user manual), etc.



Figure 1



Figure 2



Figure 3



Figure 4

- 3. Start Instrument:** Connect one end of the power adapter to the power cord and the other end to the power port of the instrument, as shown in Figure 3. Once the instrument's display screen lights up and enters the software interface, as shown in Figure 4, you can start using the instrument. For detailed instructions, please refer to the user manual.

4. Installation and Troubleshooting:

- If the instrument screen doesn't light up after connecting the power
 - i. Check if the power cord is securely connected
 - ii. Verify that the connected power source matches the requirements stated on the nameplate.